**HAYER & BOECKER**

**DIE DRAHTWEBER**

## **TYLER Ro-Tap<sup>®</sup> RX-29 / RX-30**

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### RO-TAP Repair Parts Kit

Part No.	Artikelnummer	Description
10121	562454	Minor Maintenance Kit (RX-29 & RX-94)
10122	562461	Major Repair Kit (RX-29) <i>includes #10121 kit</i>
10123		Major Repair Kit (RX-94) <i>includes #10121 kit</i>

Bild 1.1


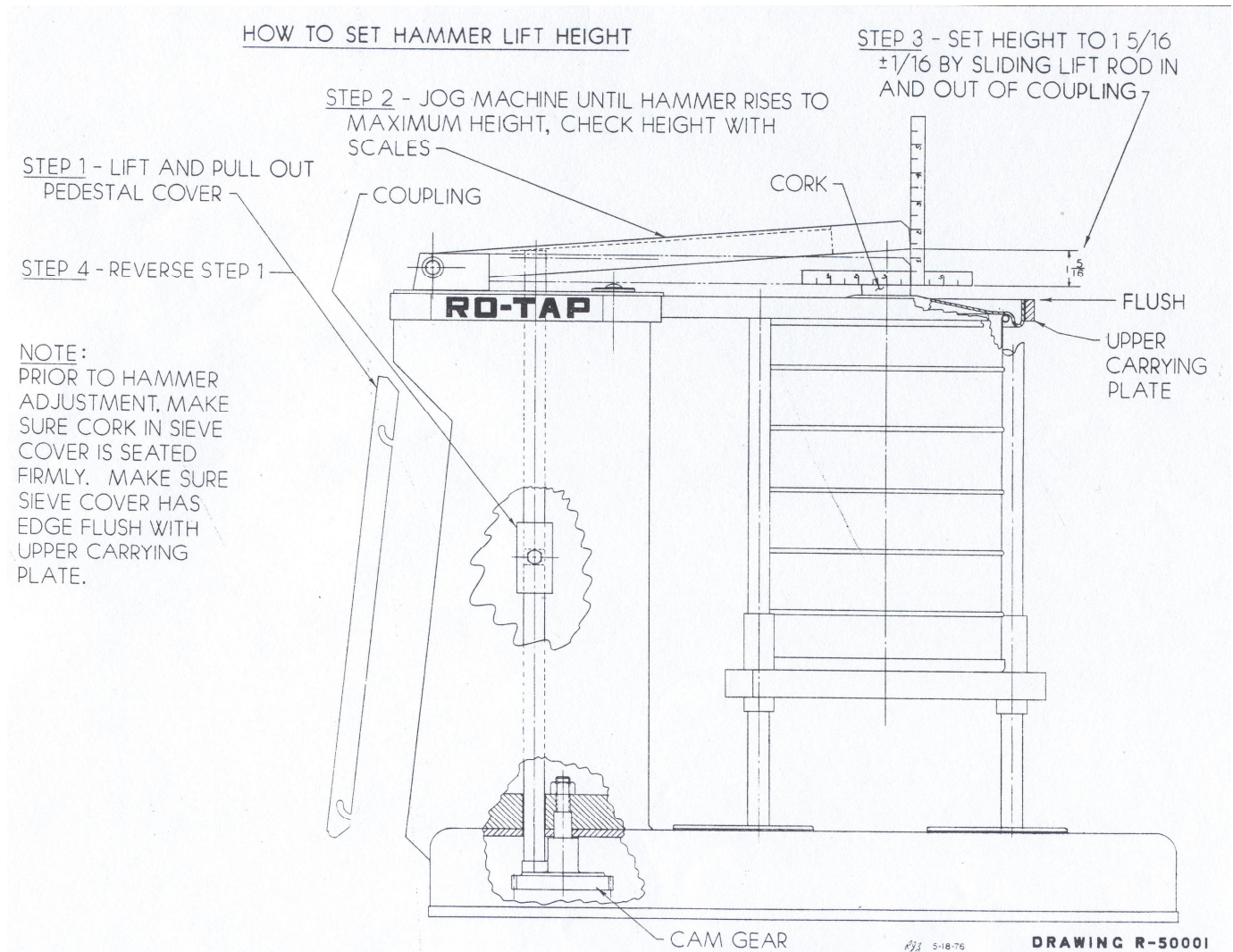
Bild 2.1
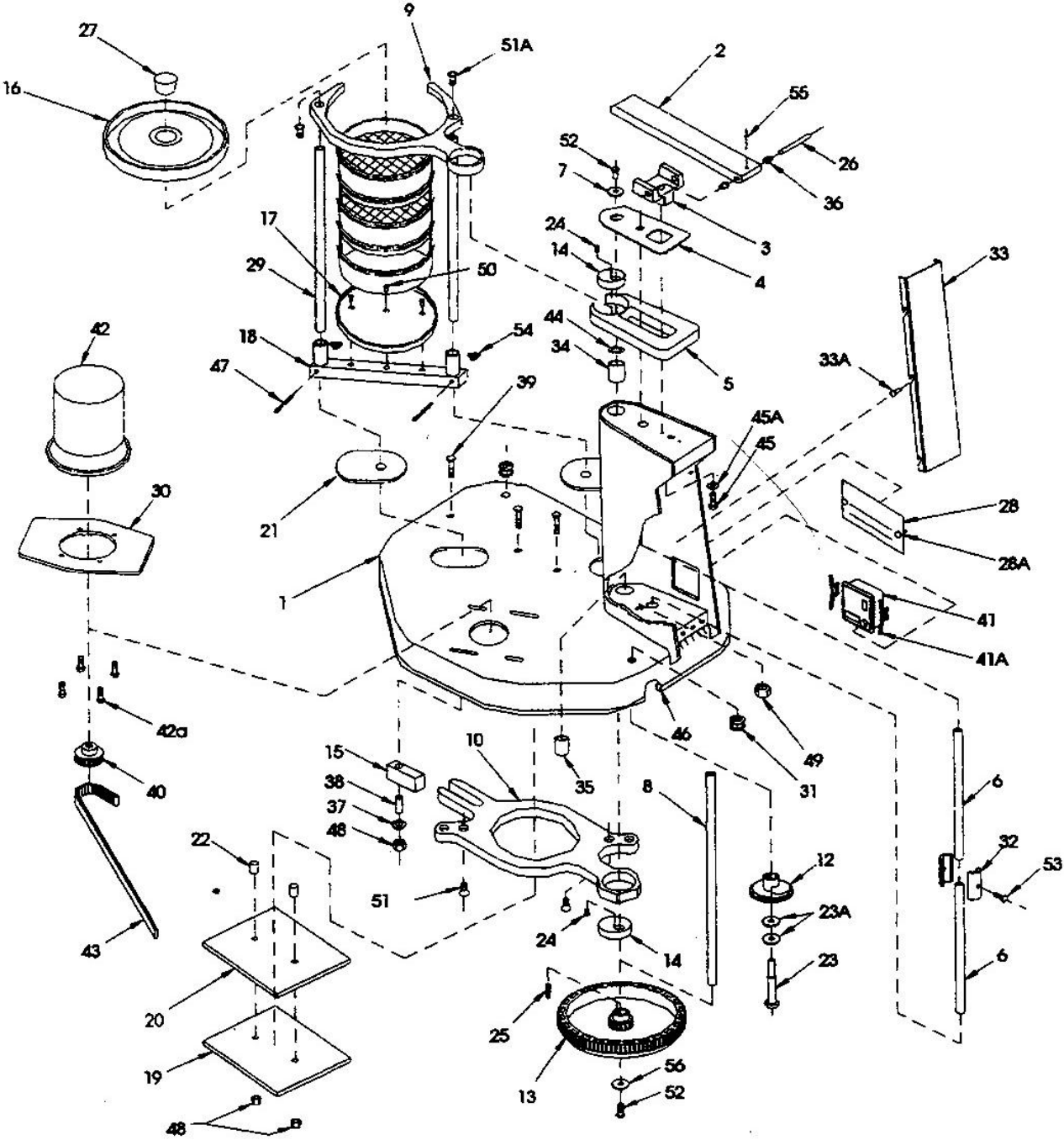
## Index

		Page
	Technical data / manufacturers proof	17
1.	Introduction	18
2.	Safety	18
2.1	Danger indications and symbols	18
2.2	Application references	19
3.	Assembly	19
4.	Handling	20
5.	Setting the hammer lift height	20
6.	Starting the shaker	21
7.	Performing the sieve analysis	22
8.	Determination of bulk volume	22
9.	Test sieves	23
10.	Sieving with the Ro-Tap <sup>®</sup> Test Sieve Shaker	23
11.	Operation time of sieving analysis	24
12.	Weighing the sieve residue	24
13.	Information about the proper use of test sieves	25
14.	Maintenance	25
15.	Guarantee	25
16.	Spare parts	25
17.	List of spare parts and drawing	26-29
	Annex: EC-Attestation of conformity, Table of test sieves and accessories	

## Technical data:

Operation voltage:	230 Volt, 50 Hertz or 110 Volt, 60 Hertz with transformer
Electronic Timer:	0-99 minutes
Power consumption:	2,9 Ampere (230V, 50Hz), 4,4 Ampere (110V, 60Hz)
Revolutions / minute:	278 +/- 10
Tappings / minute:	150 +/- 10
Weight of sieving material:	max. 3 kg (depends on bulk volume)
Dimensions test sieve shaker:	710 x 530 x 635 mm (W x D x H)
Sound Absorbing Cupboard.:	890 x 610 x 740 mm (W x D x H)
Weight Test Sieve Shaker:	approx. 80 kg
Sound absorbing cupboard:	approx. 46 kg.
Emitted sound:	The emitted sound of the machine in action has been determined with a comparative sound measurement according to DIN 45635-01 K11. The equivalent continuous sound level A is: $L_{pAeq} = 70$ (dB)

**Country of origin:** Made in U.S.A.

### Manufacturer:

W.S. Tyler Incorporated  
8570 Tyler Blvd.  
44060 Ohio / U.S.A.

### Distributor in Europe:

Haver & Boecker  
Ennigerloher Straße 64  
D-59302 Oelde

Phone.: + 49 - (0)2522 - 300  
Fax.: + 49 - (0)2522 - 30404  
E-Mail: PA@haverboecker.com

## 1. Introduction

**Only install and start the test sieve shaker after having read the following operating instructions.**

The operating instructions give important information about the secure, appropriate and economic operation of the test sieve shaker. They should be respected in order to avoid any dangers as well as to maximize the reliability and durability of the machine.

The safety instructions and regulations of prevention of accidents, as well as the different protective recommendations of the technical associations and employer's liability insurance associations, have to be strictly observed when using the machine.

**All safety instructions within these operating instructions have to be followed!**

**Always keep the operating instructions with the machine.**

If used properly, the construction of the machine guarantees that it runs and can be operated without any risk to persons.

**The guarantee does not apply to damages caused by improper use due to a non-respect of the operating instructions.**

All pictures and drawings in these operating instructions are general illustrations. They do not reflect the construction in detail.

The technical data and dimensions are without obligation, claims cannot be derived.

We reserve technical changes of the machine described below, without changing the operating instructions.

All rights on these operating instructions, the added drawings and other documents remain to HAVER & BOECKER. So are the rights on disposition, copying and passing on, including trademark registering.

## 2. Safety

### 2.1 Danger indications and symbols

The following symbols in these operating instructions mark instructions for a safe operation with the machine.

**Follow closely the instruction marked by this symbol in order to avoid any impairments of the machine or other goods.**

**Handle with particular attention to avoid health hazard.**

## **2.2 Application references**

The machine has been constructed according to the actual technical standards and the recognized safety specifications. Although, there might be dangers for the user or third persons or impairments of the machine and other goods.

Only use the machine in perfect technical constitution. Follow the operating instructions conscious of safety and danger. Repair defects at once by authorized and qualified persons!

The machine has exclusively been developed for grain sieving to determine particle distribution. Any other or further application is not as agreed upon. The manufacturer/supplier will assume no responsibility for damages as a result of improper use. The risk is on the user's side.

**For wet-sieving the RO-TAP needs a special "wet-kit"! Costs and information about this additional equipment upon request.**

## **3. Assembly**

Place the sound absorbing cupboard on a concrete foundation or a heavy, wooden work bench.

Install the RO-TAP to the middle of the cupboard. The rubber lining of the test sieve shaker foot prevents the machine from gliding.

Connect the mains plug of the test sieve shaker with the socket (inside, at the back of the cupboard).

**When installing the machine please take care that it is not accessible from downside. As the RO-TAP is open at the bottom, there is a risk of injury.  
If the machine is installed as described, there is no risk of injury.**

## 4. Handling

Once the proper test sieves have been selected, they should be stacked (one on top of the other) with the largest opening sieve on the top of the stack. A bottom pan should be added under the smallest opening sieve to collect “fine” material passing through the last sieve. The complete test sieve stack is placed on the sieve support plate (17) of the test sieve shaker. During its installation the hammer should be turned aside. After the sieving material has been filled, a sieve cover with a cork piece should be placed on top of the stack.

Adjust the sieve support plate by the sieve support clamp bar (18) in such a way that the sieve cover (16) has top edge flush with the upper carrying plate (9). The sieve cover must keep this height to make sure that the hammer drops from the right height to guaranty the precision of the sieve analysis.

## 5. Setting the hammer lift height

The hammer has been pre-adjusted by the manufacturer. If the sieving result is not satisfying the hammer setting should be re-adjusted. The hammer is adjusted in such a way, that it touches the cork piece in the sieve cover (16) as shown in fig. 1.1.

**NOTE: Prior to hammer adjustment, make sure cork in sieve cover is seated firmly. Make sure sieve cover has top edge flush with upper carrying plate.**

**Warning! Before opening the machine pull out the mains plug.**

- STEP 1      Remove pedestal cover.
- STEP 2      Jog machine until hammer rises to maximum height or use a box wrench on the hex head screw, located on top of the pedestal. Check height with scales.
- STEP 3      Set height to  $1\text{-}5/16 \pm 1/16$ " by loosening screw on coupling and adjusting lift rod.
- STEP 4      Tighten screw on coupling.
- STEP 5      Replace pedestal cover.

## 6. Starting the shaker

Before connecting the mains plug with the power supply make sure that the tension of the motor corresponds to your mains voltage.

The shaker can only be started when the sound absorbing cupboard is closed.

**Pay attention to the mains voltage!**

There are three ways to set the sieving time. It can be set in seconds, minutes or hours. To do so press the Mode key until the LED Sek/Min/Std shines at the desired position.

- Seconds: Indication of seconds
- Minutes: Indication of minutes and seconds
- Hours: Indications of hours and minutes

Main switch

1. Turn the **main switch to ON**
2. The display appears black. Pressing any key activates the indication.
3. Choose the desired sieving time with the arrow key.
4. Start the sieving by pressing the Start/Resume key.
5. A permanent beep announces the end of the sieving and has to be acknowledged with the Stop key
6. Should the sieving be interrupted or stopped, press the Stop key and restart with the Start/Resume key.

The timer can also indicate **time in general**.

**Setting the clock:** Press the Set/Display key until the LED Set shines. Select the time with the arrow keys. Press the Set/Display key once again, the time is set. To read the time press the Set/Display key shortly again. The time is indicated. After a short time the time automatically returns to the indication of the sieving time.

If the main switch is set to OFF, the machine stops. To continue the sieving set the main switch to ON and press the Start-Key. By opening and closing the cupboard's door the same function is achieved. If the sieving has to be interrupted for a longer time, the sieve time has to be re-adjusted.

## **7. Performing the sieve analysis**

In order to receive precise sieving results, the kind of procedure should be determined before starting the sieve analysis. The bulk volume, the feeding quantity, the sieving time and the desired sieving precision should be pre-determined. The RO-TAP has been constructed for dry sieving. For wet sieving with the RO-TAP a special wet-sieving-kit is available.

## **8. Determination of bulk volume**

The optimum bulk volume can be determined by regarding several factors: the weight of bulk material and its material properties such as sieving capacity, particle size and distribution. For example a sieve analysis of a solid material with a coarse particle distribution needs a feeding with bulk material of about 25 to 100 grams.

Often too large feeding quantities are filled in. To receive uniform results it is better to fill in smaller, well taken sample quantities. During a reliable sieve analysis each solid particle must have the possibility to be retained or to pass the sieve openings to the smaller ones of the following sieve. In a test sieve set of six or seven sieves a small particle must repeat this operation six or seven times. If the test sieves are overloaded a small particle may not have the chance to drop down to the right sieve bottom. However the feeding quantity has to be large enough to enable the first test sieve to retain enough particles, that are representative for the good to be analyzed.

The weight of the sieving material should be limited to a size that no sieve in the sieve set shall be overloaded. Overloading of test sieves may arise when sieving analyses are effectuated with a high degree on cut-point-close particles in the sieving material. To avoid overloading the capacities of the test sieves should be taken into account, especially the capacity of the sieve that has to retain the largest residue. Any overloading leads to uncertain results, because a lot of meshes may be blocked by particles.

The following procedure is recommended to determine the feeding quantity for test sieves: Divide the material to be analyzed with a sample splitter into samples with different weights: approx. 25, 50, 100, 150 and 200 grams. Analyze these samples on the test sieve set for about 5 minutes. When comparing the results it should be obvious, which feeding quantity leads to an optimum result. If the results, defined in percent of the residue, of a 100 gram sample are the same as the 50 gram and 150 gram sample and show a smaller mass fraction on the finest test sieve, 100 gram is the right feeding quantity to perform a sieving analysis.

Cut-point-close particles are particles, that show a particle size approaching the cut point of a test sieve. To achieve correct analyzes this test sieve should be fed slightly to enable each of the cut-point-close particles to meet the openings as often as possible.

## 9. Test Sieves

Test sieves with 200 mm or 8" (203 mm) diameter can be used for most of the sieving analyses. The standard effective height is 50 mm = 2". Two test sieves with half effective height (25 mm = 1") can be used, too. If ever possible test sieves with standard effective height should be used. When disassembling test sieve sets with half height test sieves sieving material may be spilled, sieving results are not reliable.

## 10. Sieving with the RO-TAP Test Sieve Shaker

The RO-TAP Test Sieve Shaker reproduces the circular and tapping motion given to test sieves in hand sieving, but with a uniform, mechanical action.

The results of the following sieving analysis with the RO-TAP prove its superiority to hand sieving. A sample of a fine powdered low degree copper ore, cluster-free with a lot of "slime" or dust particles has been chosen and divided to a 100 gram sample with a sample splitter.

Three sieving analyses of 20 minutes on two RO-TAP Test Sieve Shakers have been executed with the same sample for each of the three analyses. Analysis 1 and 2 have been done with RO-TAP "A", analysis 3 with RO-TAP "B". The results underline that all data can be compared.

Aperture Width in Micrometer	A.S.T.M. No.	Tyler Mesh	Test No. 1 RO-TAP A Mass fraction in %	Test No. 2 RO-TAP A Mass fraction in %	Test No. 3 RO-TAP B Mass fraction in %
600	30	28	33.4	33.5	33.7
425	40	35	18.1	18.0	18.0
300	50	48	11.0	11.1	11.1
212	70	65	8.2	8.2	8.2
150	100	100	7.2	7.3	7.2
75	200	200	9.3	9.4	9.4
	sieve pan		12.8	12.5	12.4

The results of the three sieving analyses nearly are the same. Tests No. 1 and Tests No. 2 show that the machine guarantees reproducible analyses with the same sieving material. Comparing the analyses No. 1 and No. 2 with analysis No. 3 shows that the sieving action of the two machines is nearly identical. Up to six test sieves with full effective height or up to 13 test sieves with half effective height can be used in a sieve set.

When using intermediate pans three different material quantities can be analysed with full effective height test sieves and up to seven with half effective height test sieves at the same time.

## 11. Operation time of sieve analysis

The necessary time for performing a sieve analysis with the RO-TAP depends on the kind of sieve analysis. In most cases, for example for a normal product control, three to five minutes are enough to receive the required data. Materials that are difficult to sieve may require sieve times of about 10 to 30 minutes.

When performing sieve analyses to check, whether the analysed product corresponds to a certain specification or not, longer sieving times may be necessary. Everybody concerned about should agree to the same procedure in order to receive comparable and reproducible results.

To determine the sieving time it is recommended to take three or four samples from a sample splitter with a feeding quantity that has been proved before. One of these samples should be sieved for five minutes, another for ten minutes, another for 15 minutes and a fourth one for 20 minutes. After the results have been fixed in a table, the best sieving time for uniform results shall be visible. If the result of the following sieving period does not differ more than 0.5 % to 1 % the sieving is finished.

## 12. Weighing the sieve residue

After the test sieving is finished, the whole test sieve set should be weighed by the scale to survey the data. Weighing should always be in gram with a scale with a measuring range of at least 500 gram. A sensibility of 0.1 gram is recommended.

If there are several sieve pans the residues should not be thrown away as long as the weighing has not been finished. The same procedure should be done with the test sieves in the sieve set. The sieving material retained in the sieve pan by the finest test sieve bottom has to be weighed, too, in order to receive the total weight and the proportionate in percent. The sum of the different fractions is compared to the feeding weight. The total sum of the sieving material retained on the different sieve bottoms (residue) and the residue in the sieve pan should correspond to the feeding quantity.

When using a further sieve pan (collector) to weigh the residues the material remained on the coarsest test sieve should be given first into the pan. With a smooth brass or nylon brush the downside of the sieve bottom is then brushed carefully by circular motion, without damaging the wire cloth.

In almost every cases the cut-point-close particles blocked in the meshes, can be moved out of the sieve by this dry brushing method. The test sieve border then can be tapped carefully by the brush stick to loosen last remaining particles.

### 13. Information about the proper use of test sieves

The American Society for Testing and Materials has published an information about the sieve analysis (test sieving) (STP-477).

W.S. Tyler Inc., Mentor / Ohio, describes test sieving in the Bulletin No.53. Whenever you need any particular assistance in the right handling of the RO-TAP Test Sieve Shaker and test sieves, please contact the TYLER department for laboratory equipment.

### 14. Maintenance

**Warning! Before opening the machine pull out the mains plug.**

The RO-TAP Test Sieve Shaker, Model B, requires only minimal periodic lubrication. The units incorporate plastic- and oil-impregnated bronze bearing materials, which are self-lubricating. A few drops of light oil is all that is necessary to prevent drying of the parts. Application of the same lubricants will suffice if a squeak or a drag should develop in the mechanism.

### 15. Guarantee

All functions of the sieving machine are checked and adjusted before its delivery.

The machine guarantee is for **24 months**.

**Test sieves are excluded from guarantee!**

We are only liable for defects resulting from flaws or faults in material. The guarantee does not apply to damages caused by transport. Repairs resulting by improper handling or wear due to operation are excluded from guarantee. The guarantee is void, if interventions such as variations or improper repairs are effected which are outside our control.

### 16. Spare parts

A list of deliverable spare parts is attached. Original Tyler Spare Parts are supplied by

**Haver & Boecker**

**Ennigerloher Straße 64**

**D-59302 OELDE Westphalia**

**Phone: 02522-30619 GERMANY**

**+49-2522-30619 ABROAD**

**Fax: +49-2522-30404**

**E-Mail: [PA@haverboecker.com](mailto:PA@haverboecker.com)**

## 17. List of spare parts and drawings


Item	Part No.	Article-No.	Description
1	107770	565004	Base
2	R-20029	565011	Hammer
3	R-20023	565028	Hammer Block
4	R-30019	565035	Sheet Guard
5	R-30010	565042	Rotating Guard Assembly
6	R-10036	565059	Lift Rod
7	R-10038	565066	Upper Main Shaft Washer
8	R-20027	565073	Main Shaft
9	R-30008	565080	Upper Carrying Plate (plastic)
10	R-30009	565097	Lower Carrying Plate
12	R-20039	565105	Cam Gear
13	R-30015	565112	Timing Belt Pulley
14	R-10034	565129	Eccentric Disc
15	R-10032	565136	Block
16	R-30007	565143	Sieve Cover
17	R-30006	565150	Sieve Support Plate
18	R-30023	565167	Sieve Support Clamp Bar
19	R-20019	565174	Backup Plate
20	R-20020	565181	Bearing Plate
21	R-20033	565198	Shield
22	R-10028	565206	Tube Spacer
23	R-10029	565213	Cam Shoulder Screw
23A	106582		1/2" I.D. Shim
24	R-10042	565220	Main Shaft Key to Eccentric
25	R-10039	565237	Main Shaft Key - Lower
26	R-10035	565244	Hammer Pin
27	R-10066-A	565251	Cork Plug - Material Cork
	R-10066-B	565268	Cork Plug (optional) Material Rubber
28	108184	565275	Name Plate
29	R-10033	565282	Tie Rod
30	R-30018R	565299	Motor Adaptor
31	ZA11167	565307	Grommet (replaces P-1549)
32	R-10030	565314	Lift Rod Coupling (2 1/2" lg.)
33	R-40011	565321	Pedestal Cover
34	R-10061	565338	Flange Bearing
35	R-10062	565345	Bearing
36	R-10063	565352	Flange Bearing
37	ZF-10168	565369	Thrust Washer (repl. R-10065)
38	R-10065	565376	Steel Bushing
39	ZF-10174	565383	3/8-16 NC x 2.25" Carriage Bolt (replaces P-1555-B6 Bolt)
40	R-10070	565390	Sprocket-14 teeth (1800 RPM) w/Set Screw
	R-10048	565408	Sprocket-17 teeth (1500 RPM) w/Set Screw
	R-10049	565415	Sprocket-21 teeth (1200RPM) w/Set Screw
41	R-40037	560175	Electronic Timer 50 Hz
41A	R-40040	562555	Timer Mounting Clip (2 are requested)
42	P-1500-23	565453	Motor-Standard 115/230V 60/50Hz 1 Phase
42A	ZF-10183		3/8-16 x 1.00" long Hex Head Screw
42B	ZF-10168		3/8 I.D.Flat Washer
43	R-10058	565460	Timing Belt

44	R-10055	565477	Shim
45	ZZ-10068	565484	1/2-13x2.50" long Hex Hd. Bolt (repl. P-1554-C10)
45A	ZZ-10323		1/2 I.D. Lock Washer
46	R-10052	565491	Rubber Channel (7"long)
47	R-10079	565509	Clamp Screw
48	ZF-10221	565516	3/8" Lock Nut (replaces P-1548-2)
49	ZF-10222	565523	1/2" Lock Nut (replaces P-1548-1)
50	ZF-10231	565530	10-32x"50" long Flat Head Socket Cap Screw (replaces P-1573-2)
51	ZF-10241	565547	3/8-16x .75" long Flat Head Socket Cap Screw (replaces P-1573)
52	114162	565554	3/8-16x .62" long Hex Hd. Screw
53	ZF-10251	565561	5/16-24x1.25" long Socket Head Cap Screw (replaces P-1573-4)
54	ZF-10261-B	565578	3/8 - 16 Wing Nut (replaces P-1526-B)
55	ZA-10148	565585	#10-24x.25" Socket Hd. Set Screw
56	ZF-10271	565592	3/8 I.D.x 1.50" O.D. Fender Washer
60	R-10018		Timer Cord w/Plug
61	R-10019		Cable Tie Mounting
62	R-10112		Straight Connector (APC-050)
66	107764		Flexible Conduit
67	R-10076		Wire Clip
68	ZZ-10011		Screw for Wire Clip & Ground Screw (repl. P-1569-A)


### RO-TAP Repair Parts Kit

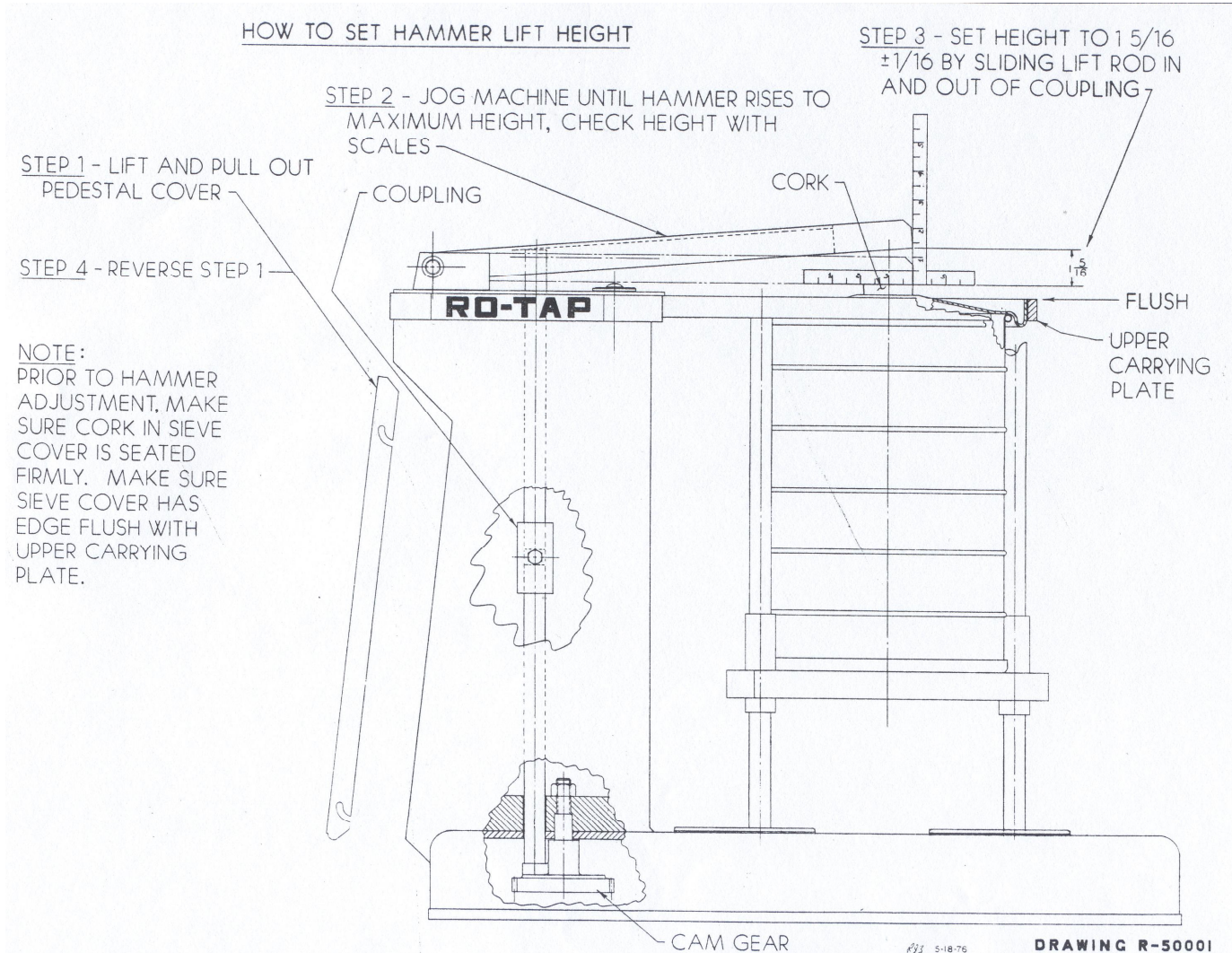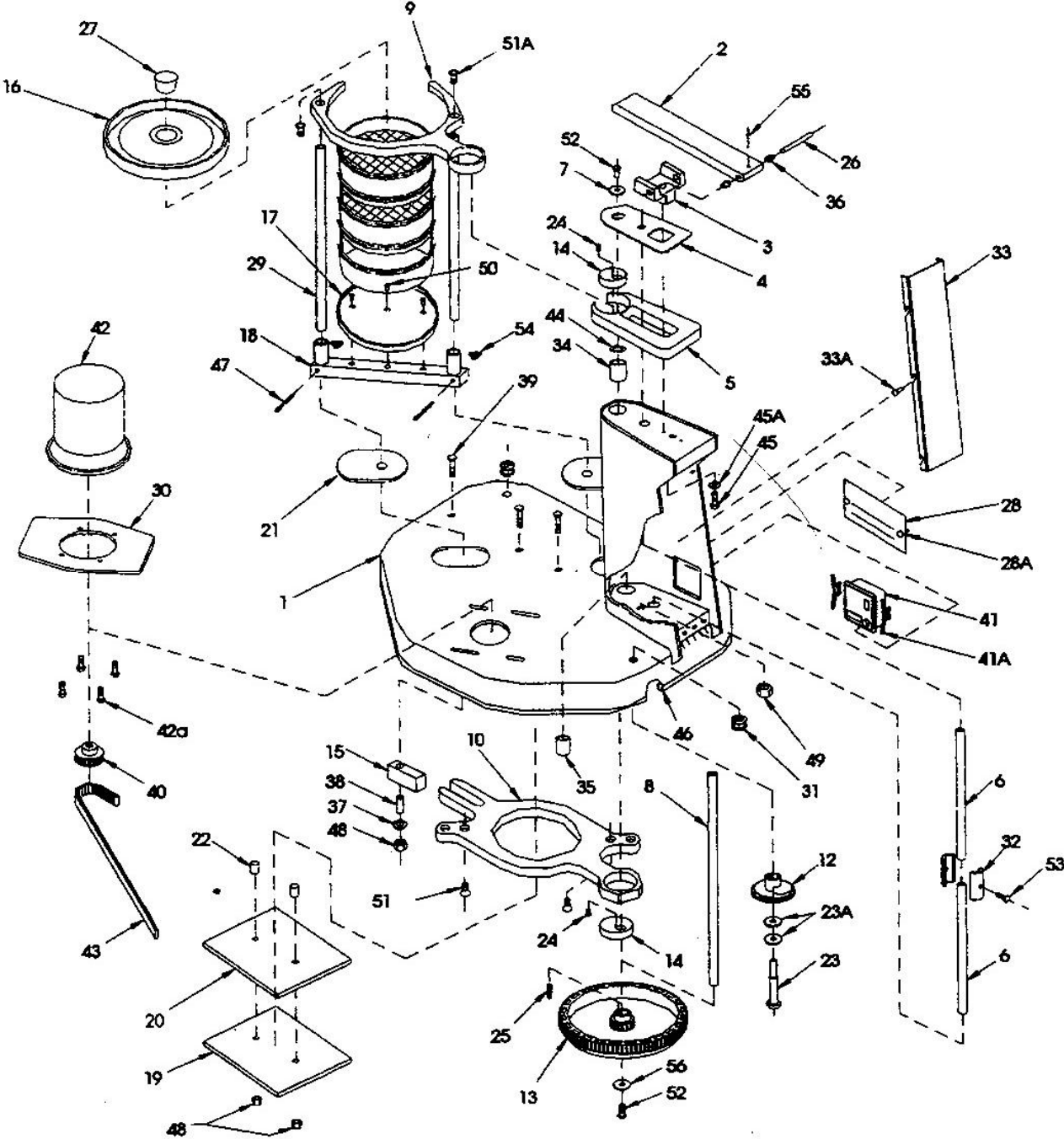Part No.	Article- No.	Description
10121	562454	Minor Maintenance Kit (RX-29 & RX-94)
10122	562461	Major Repair Kit (RX-29) <i>includes #10121 kit</i>
10123		Major Repair Kit (RX-94) <i>includes #10121 kit</i>

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